

Cytocam - Installation

1. Please check your operating system to see if you are using 32 bit or 64 bit system. To do this:

Press start (or the Windows icon)



Then go to Control Panel > System and Security > System

Under System type you will see 32 bit or 64 bit operating system.

2. Run the appropriate (32 bit or 64 bit) Camera driver application: - uEyeBatchInstall:

driver files	05/05/2015 11:16	File folder	
GigE Firmware	05/05/2015 11:16	File folder	
uEyeBatchInstall	24/10/2014 11:50	Application	216 KB

3. A Command box will appear. Press 2 (for USB drivers) then Enter.

Wait a few seconds for confirmation of driver installation. Press Enter to continue:

```
C:\Users\Magnar\Documents\Micropix\IDS\Camera drivers\64 Bit\uEye_BatchIn...
Please choose one of the following options or start the program with
the command line switch in the correspondent bracket.
The program will also copy the IDS Camera Manager to %WINDIR%\system32.
This tool can be started from the windows control panel.
1 - Install the uEye USB and Gigabit Ethernet drivers      (</all>)
2 - Install the uEye USB driver                          (</usb>)
3 - Uninstall all uEye files                             (</uinst>)
4 - Uninstall all uEye files and clean registry          (</clean>)
   ! We point out that you use this option on your own risk !
5 - Show customized command line switches
6 - Exit
Select: 2
uEye USB driver installed correctly
Press enter to continue
```

4. Run the CytocamInstaller32 application in the Cytocam folder - follow the instructions:

CytocamInstaller32	27/04/2015 16:50	Windows Installer ...	8,046 KB
setup	27/04/2015 16:50	Application	419 KB


5. Restart your PC / laptop to complete the installation.

6. Plug in your camera and wait a few seconds for the drivers to load.

Please see next page for licensing instructions.

7. Start Cytocam.

You will be prompted to load a license file (if not, click on the Micropix logo in the top left of the screen).

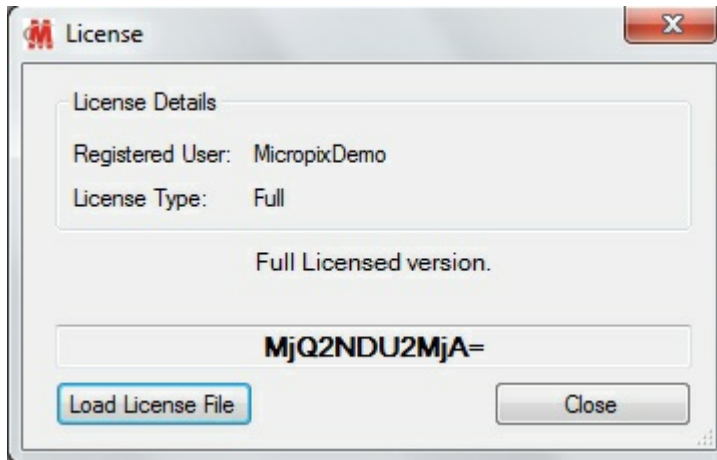


1. Click load license file.

2. Navigate to your license file (it is supplied on a memory stick with your camera). Please email us the Code if you misplace your license file.

3. Click OK.

4. You should get a success pop-up:

5. Click OK

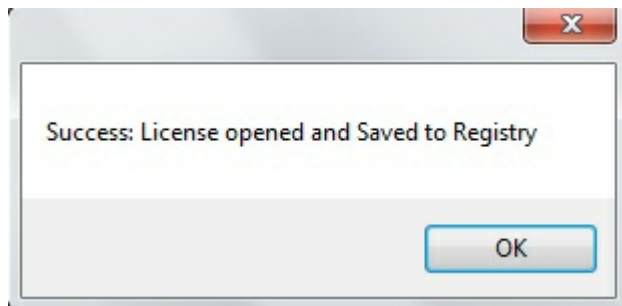

8. Installation is complete.